


HAMAMATSU

PRELIMINARY DATA
MAY 1998

MCP ASSEMBLY WITH CENTER HOLE FOR REFLECTRON MS F4294-09

Ideally suited for reflectron mass spectrometer

The F4294-09 is an MCP assembly developed for reflectron mass spectrometers. In spite of having a center hole for beam line, the F4294-09 eliminates ringing phenomena which have been a problem with conventional MCP assemblies with a center hole. The F4294-09 thus makes high-speed ion detection possible. This allows a significant size reduction in the TOF mass spectrometer and a dramatic improvement in mass resolution.

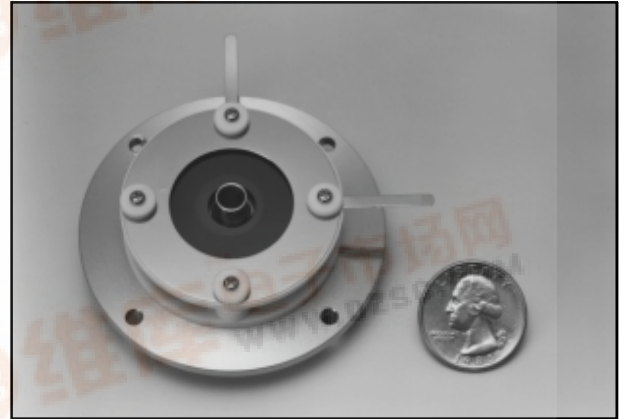

TMCPF0090

FEATURES

- Center hole for beam line
- High-speed response
- Good output waveform with negligible ringing

APPLICATION

- Reflectron mass spectrometer

SPECIFICATIONS

GENERAL

Parameter	Value	Unit
Assembly Outer Diameter	69.0	mm
Effective Diameter	27.0	mm
Maximum Height	26.4	mm
MCP Channel Diameter	12	μm
Number of MCP Stages	2	-
Center Hole Diameter	6.0	mm
Center Dead Space Diameter	12.0	mm

MAXIMUM RATINGS (Absolute values)


MCP Supply Voltage (In-Out)	2	kV
MCP-Out to Anode Voltage	0.5	kV
Electric Potential at Each Terminal	MCP-In: -2.5 MCP-Out: -0.5 Anode: GND	kV
Operating Vacuum Condition	6.7×10^{-4} (5×10^{-6}) Max.	Pa (Torr)
Baking Temperature	- *	°C
Baking Time	- *	h
Baking Vacuum Condition	- *	Pa (Torr)

* Baking cannot be performed.

CHARACTERISTICS (at 1.3×10^{-4} Pa (1×10^{-6} Torr), $T_a = +25^\circ\text{C}$)

Gain at 2.0kV	1×10^6	-
Plate Resistance per MCP	10 to 100	MΩ
Dark Count at 2.0kV	5	cps/cm ²


Figure 1: Typical Output Waveform

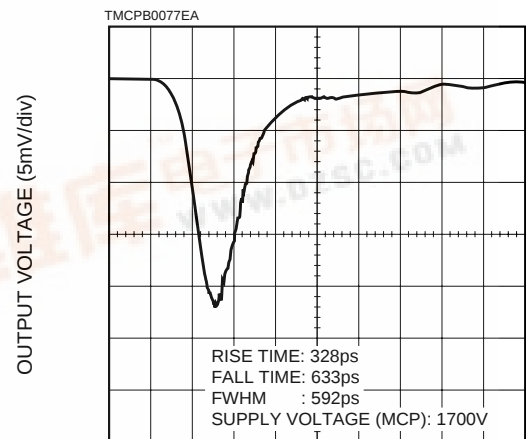

RESPONSE TIME (0.5ns/div)

MCP ASSEMBLY WITH CENTER HOLE FOR REFLECTRON MS F4294-09

Figure 2: Dimensional Outlines (Unit: mm)

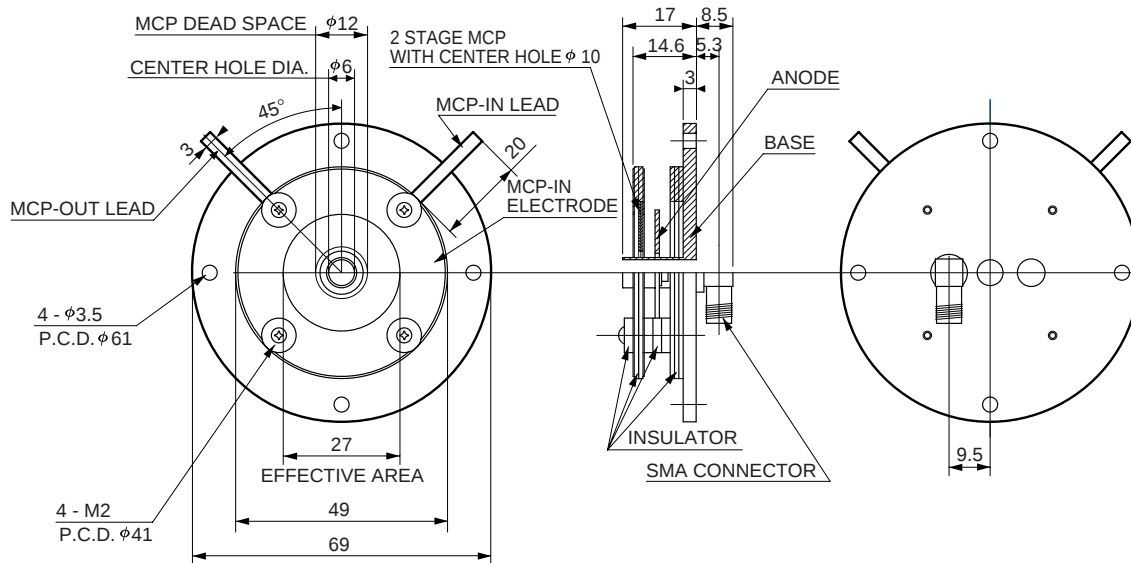

TMCPA0042EA

HAMAMATSU

HAMAMATSU PHOTONICS K.K., Electron Tube Center

314-5, Shimokanzo, Toyooka-village, Iwata-gun, Shizuoka-ken, 438-0193 Japan, Telephone: (81)539/62-5248, Fax: (81)539/62-2205

U.S.A.: Hamamatsu Corporation: 360 Foothill Road, P.O.Box 6910, Bridgewater, N.J. 08807-0910, U.S.A., Telephone:(1)908-231-0960, Fax:(1)908-231-1218

Germany: Hamamatsu Photonics Deutschland GmbH: Arzbergerstr. 10, D-82211 Herrsching am Ammersee, Germany, Telephone:(49)8152-375-0, Fax:(49)8152-2658

France: Hamamatsu Photonics France: S.A.R.L.: 8, Rue du Saule Trapu, Parc du Moulin de Massy, 91882 Massy Cedex, France, Telephone:(33)1 69 53 71 00, Fax:(33)1 69 53 71 10

United Kingdom: Hamamatsu Photonics UK Limited: Lough Point, 2 Gladbeck Way, Windmill Hill, Enfield, Middlesex EN2 7JA, United Kingdom, Telephone:(44)181-367-3560, Fax:(44)181-367-6384

North Europe: Hamamatsu Photonics Norden AB: E588natan 7, S-164 40 Kista, Sweden, Telephone:(46)8 702 20 50, Fax:(46)8 750 59 95

TMCP1010E01